

Certain hotkeys have been made available within Review to help facilitate quicker execution of certain functions from the keyboard and/or mouse without having to click additional icons or access menu options. Please note that they hotkeys are currently available only within Ponemah Review and cannot yet be augmented by the user. Please note: these hotkeys are only compatible with Ponemah v5.20.

Key Combination	Description of Action
Zoom/Graphical Navigation	
<Alt>+<S>	Synchronize Primary, Trend, and Scatter graphs with associated spreadsheet
<+> (on keypad)	Expand graph
<-> (on keypad)	Compress graph
<Ctrl>+<+> (on keypad)	Expand graph by factor
<Ctrl>+<-> (on keypad)	Compress graph by factor
<Shift>+Mouse wheel	Expand/Compress
<Ctrl>+Mouse wheel	Expand/Compress graph by a factor
Page Up	Scrolls waveform window to the left (increase in time) by screen time
Page Down	Scrolls waveform window to the right (decrease in time) by screen time
<Shift>+<F9>	Auto scale waveform in a window.
<Ctrl>+<F9>	Auto scale all waveforms in window.
<Ctrl>+<1> - <Ctrl>+<5>	Change time scale to predefined entries defined within Advanced settings of Options Application Configuration: Review.Graph.Times n
<Shift>+<6> - <Shift>+<9>	Scale one Y axis to one of the 4 defined entries (see above): Review.Graphs.High.Scalen Review.Graphs.Low.Scalen
<Ctrl>+<6> - <Ctrl>+<9>	Scale all Y axis to one of the 4 defined entries (see above)
<Ctrl>+<Z>	Undo zoom on waveform (only one level)
Mouse Wheel	Scroll left/right
<↑>	Move up one trace area – activate trace area
<↓>	Move cursor down one trace area – activate trace area
<←>	Move cursor left
<→>	Move cursor right
<Ctrl>+<←>	Move cursor the left by a factor
<Ctrl>+<→>	Move cursor the right by a factor

Key Combination	Description of Action
Add Events	
<Shift>+<A> - <Shift>+<J>	Add Event to currently selected Group
<Shift>+<Ctrl>+<A> - <Shift>+<Ctrl>+<J>	Add Event to all Groups
 on Event	Delete selected Event
Mark Placement/Positioning	
<Ctrl>+<P>	Insert Pstart and Pend Mark
<Ctrl>+<Q>	Insert QRS Marks
<Ctrl>+<T>	Insert Tend and Tpeak Mark
<Tab>	Move to next validation mark
<Shift>+<Tab>	Move to previous validation mark
<Shift>+<X>	Add Bad Data Mark
<Shift>+<F2>	Move to next Bad Data Mark
<Shift>+<F3>	Move to previous Bad Data Mark
<Shift>+<F4>	Find next change (only if Diff Marks are on)
<Shift>+<F5>	Find previous change (only if Diff Marks are on)
<Ctrl>+ <←> or <→> on Validation Mark or Bad Data Mark	Move selected mark left of right
 on Validation Mark	Delete selected mark Note: If on main Cycle mark (e.g. R-mark) deletes complete cycle
 on Bad Data Mark	Delete selected Bad Data Mark
ECGPRO	
<Ctrl>+<N>	Add New Template Cycle and analyze
<Shift>+<U>	Delete Unmatched Cycles on a Channel (User will be prompt for confirmation)