




Electrocardiogram

The Electrocardiogram Analysis Module (ECG) computes physiologically meaningful parameters on a beat-to-beat basis, from digitized multiple lead ECG data. The Multiple Lead Electrocardiogram parameters provide additional information related to QT Interval Prolongation and dispersion. The analysis functions by applying a series of logical tests to the digitized ECG signal(s), utilizing criteria selected by the user. The ECG signal can have positive, negative, or bi-phasic T waves, P waves, and QRS

complexes. The graph below represents a typical canine ECG recording and its first derivative as they would appear on the monitor. Automated validation marks for Q, R, S, End of T and Beginning of P are shown (other validation marks are also available). The validation marks provide visual, on-line verification of the accuracy of the system. The following list describes the parameters calculated by the analysis module either in real-time or during subsequent analysis.

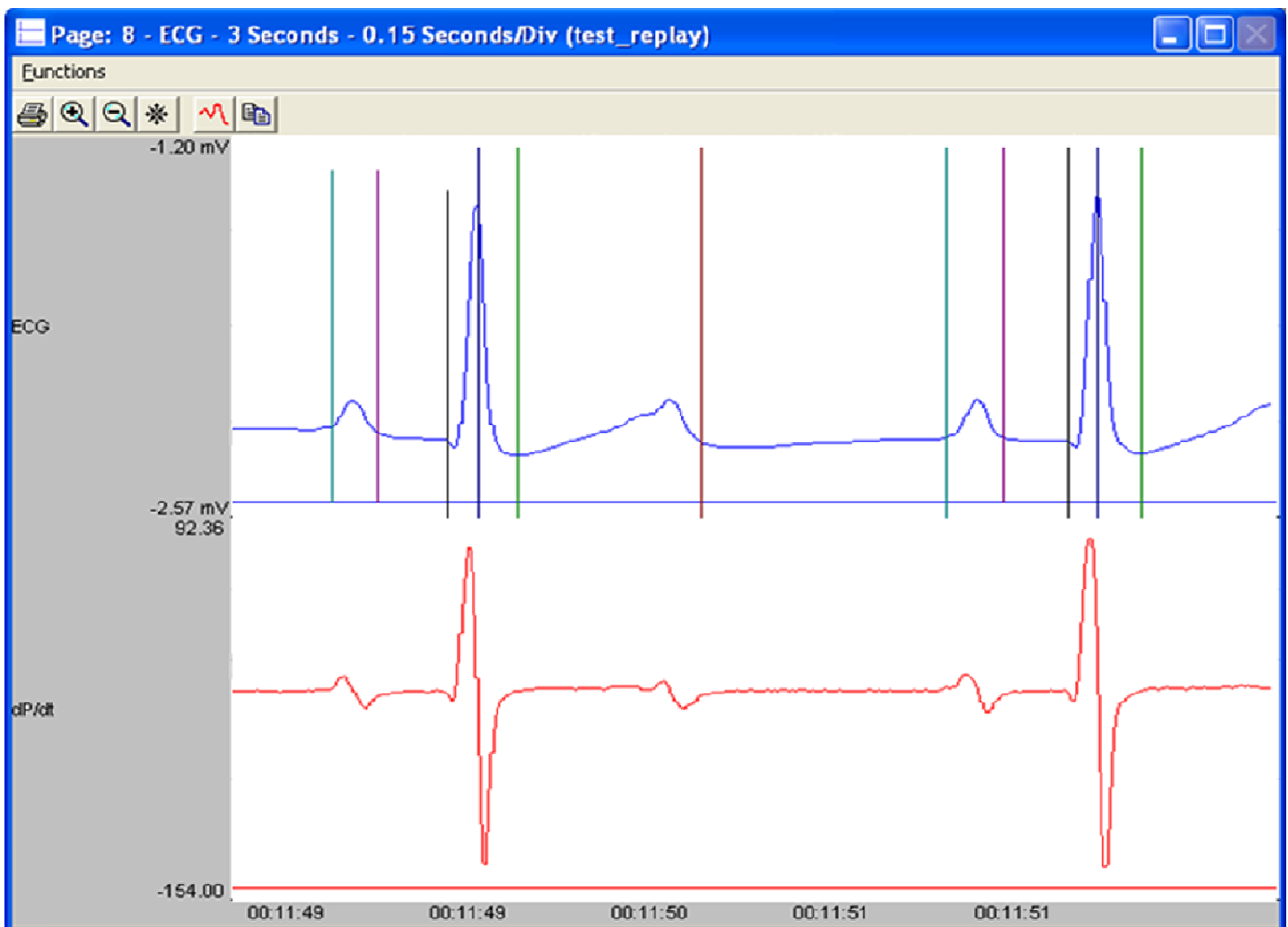

Technical Data Sheet

Model PNM-ECG100W

Electrocardiogram Analysis Module

Name	Definition
Num	The number of the cardiac cycle.
HR	The heart rate is computed in beats-per-minute.
R-H , P-H, T-H	Height of the waves from the Iso-electric level, in millivolts.
T-HN	Lowest point between the end of the S wave and the end of the T.
ST-I	Time interval in milliseconds from the S wave to end of the following T wave.
ST-E	The ST elevation, measured in "ST Measure" milliseconds after the S wave, from the Iso-electric level.
QRS	Time interval of the QRS complex.
PR-I , QT-I, RR-I	Common Interval measurements.
QAT	Q Alpha T is the time interval from the Q wave to the peak of the following T wave.
QTcb , QTcf, QTcv, and QTcm	Multiple corrected QT intervals (Bazett, Fridericia, Van de Water and Matsunaga).
EQTS, EQTSc, EQTM, EQTMcs, EQTMce, QTMc	Various cross channel calculations.
QTD	QT Dispersion.
QR-I, QRSA	QR interval, QR amplitude.
MxdV	Maximum derivative of the R wave.
T-A	Area of the T wave from the Iso-electric level.
PCt , TCt,	The number of valid waves encountered in the logging period.
QTCt	QT count.
BAD	The number of arrhythmic beats detected during a specified logging period.
GW, TW	The Good Wave counts and number (Total Wave) of good and bad complexes.
QATN	Time, in milliseconds, between the Q wave and the lowest point between the end of S and the end of T wave.
PWdth (Pwidth)	Reports the time, in milliseconds, between the start and end of the P wave.
Tpe-I	The time in milliseconds between the peak of the T wave and the end of the T wave.
T-P	The signal value at the peak of the T wave relative to the Iso-electric level.
Match, Pmatch, Qmatch, Smatch, Tmatch	Used specifically with Template Analysis.
Noise	This parameter reports an approximation of the noise level in the ECG cycle.

DSI products are not intended for the purposes of diagnosis of disease or other conditions, or in the cure, mitigation, treatment, or prevention of disease, or used as a life support device. Use of DSI products are solely for the purposes of conducting life science research.