

Installation Instructions

Ponemah Version 6.41

Prior to installing, it is recommended to review the Ponemah v6.41 Release Notes.

Ponemah v6.41 is compatible with 64 bit versions of Windows 7 and Windows 10.

NOTE:

- Installation must occur from a Windows® Administrator account.
- If using PhysioTel Digital, CLC firmware 1.29.206 is required.

To access the Ponemah install:

1. **Download it from internet:**
 - a. Go to www.datasci.com
 - b. Select the **Resources** menu.
 - c. Select **Software and Firmware Updates**.
 - d. Extract the files by **right-clicking** the .zip file and choosing **Extract**.
 - e. Browse to the extracted files and **double-click** on the **install.exe** file to display the installation options.
2. **Request a physical copy of the install media:**

The website is the primary means for delivering Ponemah. If a physical copy of the install is desired, one can be requested by contacting DSI Client Services (CSservice@datasci.com).

 - a. Place the USB drive into one of the computer's USB ports.
 - b. Navigate to the **install.exe** file and **double-click** it to display the installation options.
3. **License File Access:**
 - a. License files can be downloaded from the internet at the following FTP site:
<ftp://207.67.9.225>
User Name: client\$user
Password: ponemah\$1
 - b. Click the directory range that includes the license file serial number
 - c. Click the license file serial number to download the license
 - d. License files can also be obtained by contacting DSI Technical Support (Support@datasci.com)

To install Ponemah v6.41:

1. Click **Install Ponemah**.
2. Follow the onscreen instructions.
3. Click **Finish**.

NOTE: Previous versions of Ponemah v6.x do not need to be uninstalled.